



UC Davis MMPC-Live Protocol

Triglycerides

Version: 1.0
 Revision Date: 9/17/2012
 Replaces version: None
 Edited by: Peter Havel - UC Davis Metabolism & Metabolic Health Core

[Summary](#)
[Reagents and Materials](#)
[Protocol](#)
[Reagent Preparation](#)

Summary:

Triglycerides are enzymatically hydrolyzed by lipase to free fatty acids and glycerol. The glycerol is phosphorylated by adenosine triphosphate (ATP) with glycerol kinase (GK) to produce glycerol-3-phosphate and adenosine diphosphate. Glycerol-3-phosphate is oxidized by dihydroxyacetone phosphate (DAP) by glycerolphosphate oxidase producing hydrogen peroxide (H₂O₂). In a Trinder5 type color reaction catalyzed by peroxidase, the H₂O₂ reacts with 4-aminoantipyrine (4-AAP) and 3,5-dichloro-2-hydroxybenzene sulfonate (DHBS) to produce a red colored dye. The absorbance of this dye is proportional to the concentration of triglycerides present in the sample.

Reagents and Materials:

| <i>Reagent/Material</i> | <i>Vendor</i> | <i>Stock Number</i> |
|-------------------------|--------------------|---------------------|
| Calibrator | Fisher Diagnostics | TR43002 |
| Reagents | Fisher Diagnostics | TR22203 |
| PBS | | |
| Microplate | | |
| Platereader | | |

Protocol:

1. Reconstitute powdered reagent with only 25 ml of distilled water to make a 2X solution.
2. Add 3 μ l of calibrator and sample to each well.
3. Add 150 μ l of PBS to each well. Read at 540 nm.

IMPORTANT: Make sure not to add any bubbles to the wells when dispensing reagents, this will interfere with reading in the platereader.

4. Add 150 μ l of 2X reagent to each well. Incubate at 37°C for 5 minutes. Read at 540 nm.
5. Subtract blank readings from final readings. The assay will be linear so the unknown samples can be calculated as (Sample Absorbance \div Calibrator Absorbance) \times Calibrator Concentration.

Reagent Preparation:

PBS – ready to use

Reagent – reconstitute with distilled water to make a 2X solution